

AN AXION™ EYEWASH...



...FOR THE OFFICE,
HOME AND LAB

AXION™ eyePOD™

The AXION™ eyePOD™ is a faucet-mounted eyewash that provides professional business facilities, such as doctor and dentist offices, laboratories, and educational facilities, with AXION MSR™ eyewash capabilities in an attractive, low profile, low cost design. With a flick of the wrist your standard faucet becomes a thermally controlled AXION™ Medically Superior Response™ eyewash, and with another flick of the wrist it resets as a faucet. Now how is that for convenience?

axion™
EYEPOD™


AXION MSR™

MEDICALLY SUPERIOR RESPONSE™

A medical problem ... an elegant solution

AXION MSR™ eyewash products use an exclusive, inverted flow pattern that is consistent with first aid protocols used by medical professionals. The eyewash flows mitigate the risk of inhalation or ingestion of hazardous contaminants during eye irrigation.

The medical profession teaches and practices irrigating eyes by introducing flushing fluid at the inner corner of the eye – adjacent to the nose – letting it run across the eye to the outer edge. This is specifically done to avoid driving contaminants into the lacrimal punctum and the nasal cavity. The accepted medical practice is counter to the way that traditional eyewashes operate and can result in further injury to the splash or blown-particulate injured victim. AXION MSR™ is the only eyewash product that duplicates the methodology used by and recommended by healthcare professionals.

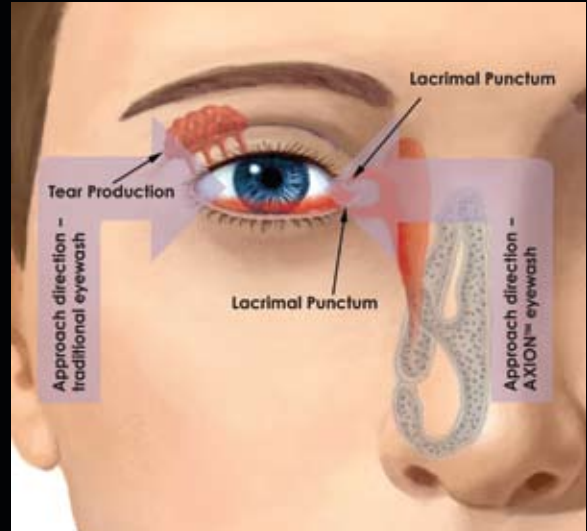


The newest addition to the AXION™ line of products, the AXION™ eyePOD™, is a faucet-mounted eyewash that provides professional business facilities with AXION MSR™ eyewash irrigation capabilities in an attractive, low cost, compact package.

The sleek, smooth, stainless steel exterior of the eyePOD™ blends in with any décor and is easy to clean and use.

Installation is a snap, with a variety of industry standard adapters included that allow the eyePOD™ to be used with standard faucets. Simply remove the existing aerator from the bottom of your faucet by unscrewing it and replace it with the eyePOD™.

The elegant eyePOD™ has been designed to withstand a lifetime of use and is certified to NSF61, ANSI/ISEA Z358.1.

As an added safety precaution, the eyePOD™ features a built-in thermostatic safety valve that will automatically shut off eyewash flows at 100°F (37°C), diverting all water out through the aerator at the bottom of the unit. This protects users from accidentally irrigating with water that is too hot. When water temperatures fall to an acceptable degree, eyewash flows are automatically restored.

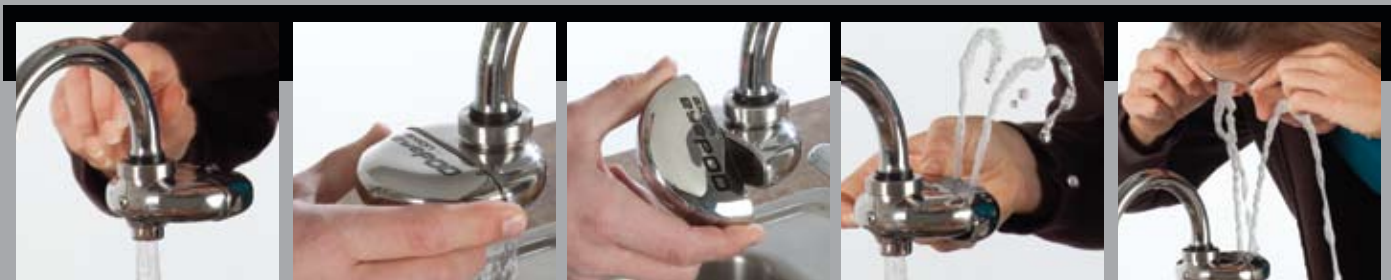

With the attractive logo facing upward, the AXION™ eyePOD™ will allow your faucet to function normally. Water passes through the eyePOD™ and exits through the built-in aerator under the unit.

Rotating the front of the eyePOD™ one-half turn (in either direction) brings the eyewash spouts to the top and automatically resets the unit to operate as an eyewash. Water is discharged from the twin AXION™ eyePOD™ eyewash nozzles for a soothing irrigation

experience. The downward orientation of the eyewash nozzles when not in use keeps them free of dust and debris accumulation... and, their antimicrobial treatment minimizes the risk of damage due to mold and mildew growth.





www.axioneyepod.com

US/Canada

Haws HQ • www.hawSCO.com • 1.888.640.4297

Mexico/Central America

Haws HQ • www.hawSCO.com • +1.775.359.4712

South America

Avlis-Haws Brazil • www.avlisvalvulas.com.br • +55 (11) 4024.2977

Asia Pacific

Haws Singapore • www.haws.sg • +65.6367.0359

Europe/Middle East/Africa

Haws Switzerland • www.haws.ch • +41 (0) 34.420.6000